Chapter 10 • Chemical Quantities

**EXPERIMENT** 



# EMPIRICAL FORMULA DETERMINATION

# PURPOSE

To determine the empirical formula of magnesium oxide.

Text Reference Section 10.3

#### BACKGROUND

Carbon dioxide (CO<sub>2</sub>), water (H<sub>2</sub>O), and ammonia (NH<sub>3</sub>) are three out of many chemical compounds that you are familiar with. Have you ever seen a compound with a formula such as Na<sub>2.3</sub>Cl<sub>3.9</sub>? In fact, such a formula is impossible. Only whole atoms, not fractions of atoms, react with each other to form products. Also, although elements may react in different proportions to form more than one compound, the proportions of atoms in those compounds will always be a ratio of small whole numbers.

An empirical formula gives the simplest whole-number ratio of the different atoms in a compound. For example, while the molecular formula for hydrogen peroxide is  $\rm H_2O_2$ , the simplest whole-number ratio of hydrogen and oxygen atoms can be expressed as HO. Thus, the empirical formula of hydrogen peroxide is HO.

In this lab, you will experimentally determine the empirical formula of magnesium oxide, the compound formed when magnesium metal reacts with oxygen.

## **MATERIALS (PER PAIR)**

safety goggles crucible crucible lid clay triangle ring stand ring support

crucible tongs
gas burner
centigram balance
magnesium ribbon, Mg F
2 pieces of exposed film

#### **SAFETY FIRST!**

In this lab, the crucible you are using will become quite hot and could cause a severe burn if handled improperly. Observe all precautions, especially those listed below. If you see a safety icon beside a step in the Procedure, refer to the list below for its meaning.



Caution: Wear your safety goggles. (All steps.)



**Caution:** Do not look directly at burning magnesium. Do not inhale the smoke produced when the magnesium is burned. (Steps 3, 4.)



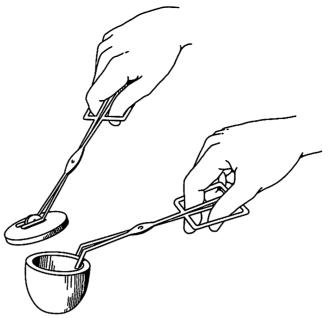


Figure 13.1



Caution: Always handle the crucible and crucible lid with crucible tongs, as shown in Figure 13.1. (Steps 3, 4.)



Note: Return or dispose of all materials according to the instructions of your teacher. (Step 7.)

#### **PROCEDURE**

As you perform the experiment, record your data in Data Table 1.



- 1. Set up the equipment, as shown in Figure 13.2. Clean a crucible and its cover with water. Dry them by heating in the hottest part of the flame for 5 minutes. Allow them to cool for at least 10 minutes. Measure and record the combined mass of the crucible and lid to the nearest 0.01 g.
- 2. Place a coiled 25-cm length of magnesium ribbon in the crucible. Measure and record the combined mass of the crucible, lid, and magnesium.







- 3. CAUTION: Do not look directly at the burning magnesium. View the reaction through the pieces of film provided by your teacher. Over a high flame, heat the uncovered crucible on the triangle until the magnesium ignites. CAUTION: Do not inhale the smoke produced. When the magnesium begins to burn, immediately cover the crucible (using tongs) and remove the burner.
- 4. After smoke production has ceased, replace the burner and continue heating the crucible. CAUTION: Do not lean over the crucible. Remove the burner and carefully lift the lid and check the reaction every 2 or 3 minutes. After heating for a total of 10 minutes, check to see if the reaction is complete. The magnesium should be wholly converted to a light gray powder, magnesium oxide. If ribbonlike material remains in the crucible, replace the burner and continue heating.

- completely (at least 10 minutes).
  - 6. Measure and record the combined mass of the crucible, crucible lid, and magnesium oxide.

5. Turn off and remove the burner. Allow the crucible to cool

- 7. Follow your teacher's instructions for proper disposal of the materials.

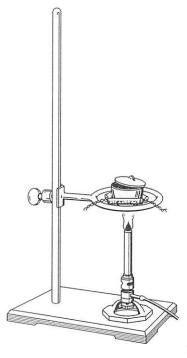


Figure 13.2

### **OBSERVATIONS**

DATA TABLE 1: MASS MEASUREMENTS				
Item	Mass			
empty crucible and lid				
crucible, lid, and Mg (before heating)				
crucible, lid, and combustion product $(Mg_xO_y)$				

#### **ANALYSES AND CONCLUSIONS**

- 1. Determine the mass of magnesium used.
- 2. Determine the number of moles of magnesium used.

**Hint:** mol Mg = (mass Mg/molar mass Mg)

- **3.** Determine the mass of magnesium oxide formed.
- 4. Determine the mass of oxygen that combined with the magnesium.
- 5. Calculate the number of moles of oxygen atoms that were used.
- 6. Calculate the ratio between moles of magnesium used and moles of oxygen used. Express this ratio in simplest whole-number form.

Name	Da	ate	Clas	SS	
11. Interpret, in terms of chemical formula for	atoms and in terms ethane $(C_2H_6)$ .	of moles, the su	bscripts in the		
			,		
GOING FURTHER	?				
Develop a Hypoth		·			
At the high temperature also be formed. Propose	at which this reaction hypothesis about t	n takes place, a he chemical nat	green-yellow soli ure of this produc	d may xt.	
	······································				
Design an Experim	ent				
Propose an experiment to your teacher's permission			e available and yo	u have	
					w-1,